



Nippon Instruments Corporation

Determination of Mercury in Virginia Tobacco Leaves (CTA-VTL-2)

Principle	Pyrolysis-gold amalgamation AAS
Related Official Methods	EPA 7473, ASTM D6722-01
Instrument Model	MA-2000
Sample (Sample Name)	Virginia Tobacco Leaves (CTA-VTL-2)

* **Data:** **n=5**

Average Value : **0.0425 mg/kg**

Reproducibility (CV%): **2.76%**

Reference (Certified) Value: **0.0480 ± 0.009mg/kg**

Remarks

- **Accurate results for Virginia Tobacco Leaves (CTA-VTL-2)**
- **Good reproducibility proved**

